

GroMed™

3 grams Soybean - Casein Digest Medium Catalog #GM3000

For the cultivation of microorganisms. Not for patient use.

Nonsterile.

Store at 2 to 30° C.

1. Combine contents of packet (approximately 3gm) with 100mL of nonbacteriostatic water to make a 3% solution.
2. Heat and stir as needed to dissolve completely.
3. Yields approximately 100mL of Soybean-Casein Digest Broth (Tryptic Soy Broth).
4. Liquid media can be inoculated with specific microorganisms or set out over night in an uncovered wide-mouth container.
5. Optional: Use in conjunction with PATT™ kits, #GM7020 or #GM7030, to perform a high-risk media fill test.

IRRITANT: Dry Powder may be irritating to eyes, skin, and respiratory system.
Avoid contact or inhalation.

This lot was found to meet or exceed the following performance criteria and applicable NCCLS (National Committee for Clinical Laboratory Standards) standards for performance.

Test Organisms		Results
<i>Escherichia coli</i>	ATCC® 25922	Growth
<i>Candida albicans</i>	ATCC® 10231	Growth
<i>Staphylococcus aureus</i>	ATCC® 25923	Growth
<i>Bacillus subtilis</i>	ATCC® 6633	Growth

Physical Characteristics

pH: 7.3±0.2 at 25°C

Fill: 3gm

Appearance: Homogeneous and free flowing

Color: Light beige

Q.I.medical, inc

Nevada City, CA 95959
(530) 265-4820 • FAX (530) 265-9416
Technical Services (800) 837-8361
www.qimedical.com

DFU020 C 05/05

