

# GroMed™ 5mL FTG in Syringes

## Fluid Thioglycollate (FTG) Growth Media Catalog #GM6105

Fluid Thioglycollate is recommended for the cultivation of microaerophilic and anaerobic microorganisms.

Store at 2° to 25°C, out of direct light, until ready to use.

1. Fluid Thioglycollate in syringes, GroMed #GM6105, are primarily intended for use with the QTMicro Systems, #TM6000 and #TM6100.
2. Follow the Directions for Use supplied with each case of QTMicros, steps #1 through #8.
3. At step #9, substitute a syringe of FTG, #GM6105, for the syringe of TSB that is supplied in the case of QTMicros.
4. At step #10, gently press the plunger on the syringe containing FTG while filling the QTMicro filter chamber. The goal is to minimize turbulence and introduction of oxygen into the media in the filter chamber.
5. At step #13, Incubation:  

USP Chapter <71> Sterility Tests: Incubate at 32.5°+/-2.5°C for not less than 14 days.
6. Follow steps #14 and #15 in the Directions for Use supplied with the QTMicro Systems.

**Q.I.medical,inc**

Nevada City, CA 95959  
(530) 265-4820 • FAX (530) 265-9416

DOC315 A 01/06